

ITEM	QUANTITY
Protective non-latex gloves (large)	2 pair
Protective shoe covers	1 pair
Protective mask with eye shield	1
Protective hair cover	1
Protective plastic apron with sleeves	1
Absorbent - 1 oz. package	1
Scraper and shovel	1
EPA-approved Sanizide Plus® disinfectant spray - 4 oz.	1
Absorbent towel	1
Antimicrobial hand-wipe	1
MSDS for absorbent	1
MSDS for disinfectant spray	1
Spill clean-up instructions	1
MSDS for towlette	1

Sharps® Biohazard Spill Clean-Up Kit


Reorder #51000 (includes mail-back pail and clean-up kit)
 Reorder #85100 (includes ship-back pail and clean-up kit)
 Reorder #50027-002 (includes 2 spill clean-up kits - no pail)



PLEASE READ COMPLETELY BEFORE PROCEEDING #51000 / #85100

Sharps® Biohazard Spill Clean-Up Kit

This Biohazard Spill Clean-Up Kit contains everything required to safely and properly clean-up biohazardous spills in accordance with OSHA and state regulations.

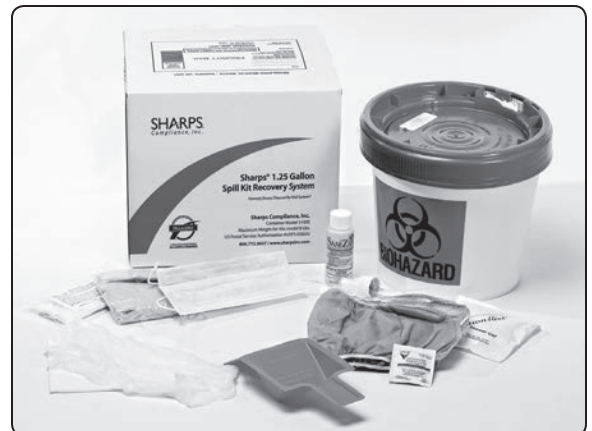

Sharps® Biohazard Spill Clean-Up Kit

Instructions for Proper Spill Clean-up

1. Secure spill to eliminate foot traffic until clean-up is complete and disinfected surfaces are dry.
2. Put on protective gloves and additional personal protective equipment (PPE) as determined necessary. (Figure 1)
3. Ensure that pail or bag is open and ready to receive waste material.
4. Open the pouch of Absorbo-Gel absorbent material and sprinkle liberally over spill. (Figure 2)
5. Allow absorbent to fully soak up the spill.
6. Use the scraper and scoop to pick up the absorbed spill material. Deposit the spill material along with the scoop and scraper into the pail or bag. (Figure 3)
7. Saturate the affected area with the SaniZide Plus® disinfectant spray and allow area to remain for time recommended on the spray bottle. Spray multiple times as necessary to completely clean the area. (Figure 4)
8. Wipe disinfected area with the paper towel. Discard paper towel in the pail or bag.
9. Remove PPE carefully and discard into pail or bag. (Figure 5)
10. Clean hands using the antimicrobial hand-wipe. Discard wipe into pail or bag.
11. Secure lid onto pail tightly or tie up bag. Wash hands with soap and water as soon as possible.
12. If applicable, follow instructions included with the #51000 or #85100 Recovery Systems to properly package and return system to Sharps Compliance for proper treatment. For more information, go to www.sharpsinc.com
13. Comply with all state and local regulations regarding proper disposal.



Figure 1

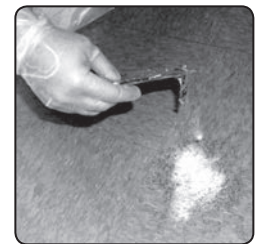


Figure 2



Figure 3

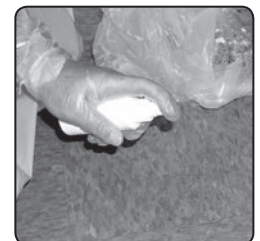


Figure 4



Figure 5

**FOR ASSISTANCE, contact Sharps at 1.800.772.5657
or www.sharpsinc.com**